How to Install SNPViewer and Visualize Data

Milcah Kigoni (m.kigoni@cgiar.org)

Sep 2018





Step 1: Download SNPviewer from the link below:

https://www.lgcgroup.com/products/genotyping-software/snpviewer/



> Genotyping software

SNPviewer is suitable for computers that are running on Windows, 2000,

Step 2: You will receive results in a zip file with three file. Unzip and get the .csv file similar to the one below . This is the SNPViewer input file

	ן \$י כי י י י י	Ŧ					000 - SNPviewer-012.015-01 - Excel										困 -	đ	\times
Fi	le Home Inse	ert Pagel	Layout Fo	rmulas	Data Review	w View	🗘 Tell me	what you wa	int to do						Kig	oni, Milcah	(ICRISAT-Nai	robi) 🖓	Share
Past	Cut Copy → Format Painter Clipboard □	Calibri B I U	 ▼ 11 ▼ 	• A A > - <u>A</u> •		N→ I III Alignment	Wrap Text Merge & Cer	Gene nter - \$ -	eral % ∮ €.c Number		itional Form atting - Tab Styles	nat as Cell ole • Styles •	Insert D	elete Format	∑ AutoSun ↓ Fill ~	n • AZY Sort & Filter • Editing	Find & Select •		^
B5 \cdot : \times \checkmark f_x 0																~			
	А	В	С	D	E	F	G	Н	I	J	К	L	М	Ν	0	Р	Q	R	
1	KBiosciences genot	typing rep	ort																
2	LGC-Genomics																		
3	Export version	1.21																	
4	Customer	Wheat																	
5	Customer number	C																	
6	Project number	C)																
7	Master plate type	96	5																
8	Order number	C)																
9	View mode	auto / ori	ginal data																
10	Viewing	Please download SNPViewer to graphically view this file or to find out more about the syntax from http://results.lgcgenomics.com/software/snpviewer/																	
11	Editing	This file maybe graphically edited and re-called using LGC KlusterCaller software from http://results.lgcgenomics.com/software/klustercaller/																	
12	Title	000 XXXC	-B01 16 pla	tes															
13																			
14	Statistics																		
15	SNP	Plate	Υ	Het	Х	NTC	Unused	Missing	Bad	Allele Y%	Allele X%	X2	Allele Y	Allele X					
16	snpTA0016	17738-01	20	2	25 44	. 2	2 5	0	(36	52	63	48	3 13	83 T		G		
17	snpTA0016	17738-02	0	2	26 66	2	2 2	0	() 14	13	85	87	2	49 T		G		
18	snpTA0016	17738-03	0	2	25 68	2	2 1	0	() 13	44	86	56	j 2	24 T		G		
19	snpTA0016	17738-04	46		0 42	2	2 5	0	(52	27	47	73	88	T 0		G		
20	snpTA0016	17738-05	38		5 46	2	2 5	0	() 45	51	54	49	69	98 T		G		
21	snpTA0016	17738-06	1	. 1	13 76	2	2 3	0	() 8	33	91	67	0	27 T		G		
22	snpTA0016	17738-07	15	1	13 63	2	2 3	0	(23	63	76	37	33	21 T		G		
	> 000 - SN	Pviewer-0	12.015-01	(+)							:	•							



Load the .csv file to SNPviewer: Visualize and export genotyping results as cluster plots



Use the guidelines here to run the to visualize the data:

https://www.lgcgroup.com/LGCGroup/media/PDFs/Products/Genotyping/Software/SNPviewer_manual.pdf?ext=.pdf